

NutriStem® hPSC XF Medium Supports Long-term Culture of Human Pluripotent Stem Cells as Clumps on Corning® Matrigel® Matrix and Single Cells on Corning PureCoat™ rLaminin-521 Cultureware

CORNING

Jeff Partridge, Himabindu Nandivada, Paula Flaherty, and Deepa Saxena
Corning Incorporated, Life Sciences, Bedford, MA USA

Introduction

Human pluripotent stem cells (hPSCs) including human embryonic stem cells (hESCs) and human induced pluripotent stem cells (hiPSCs), are typically cultured using complex media containing animal-derived components. In addition, adherent culture of hPSCs requires either a layer of feeder cells (e.g., mouse or human fibroblasts), a complex mixture of naturally derived extracellular matrices (e.g., Corning® Matrigel® matrix) or defined surfaces (e.g., recombinant Laminin). In this study, we demonstrated long-term culture of hiPSCs in NutriStem® hPSC XF medium – a serum-free, xeno-free medium – widely used for long-term growth and expansion of hPSCs. Human iPSCs were cultured as clumps on Matrigel matrix or as single cells on Corning PureCoat™ rLaminin-521 cultureware in NutriStem medium for at least ten passages. Cells exhibited typical hPSC morphology and remained undifferentiated as demonstrated by the expression of Oct-3/4 (>94%), SSEA-4 (>95%) and the absence of SSEA-1. After ten passages, pluripotency was shown by differentiation into the three germ layers and cells retained a normal karyotype.

Materials and Methods

Matrigel coated plates were prepared using Corning Matrigel hESC-qualified matrix (Cat. No. 354277). Corning PureCoat rLaminin-521 cultureware (Cat. No. 356290) was used according to manufacturer's guidelines. Human episomal iPSCs (Gibco®) were routinely maintained on Matrigel coated 6-well plates in mTeSR®1 medium (STEMCELL Technologies™).

On Matrigel coated 6-well plates, cells were clump passaged every three to five days using Cell Dissociation Buffer (Life Technologies) and seeded at a split ratio of 1:14 - 1:21 in NutriStem XF/FF medium (manufactured by Biological Industries).

Cells were seeded as single cells onto rLaminin-521 cultureware (6-well format) in NutriStem medium. Cells were passaged every four to five days using Accutase® and seeded at a density of 50,000 cells/cm².

At each passage, the morphology of the cells and the cell yields (PureCoat only) were monitored. After ten passages, cells were karyotyped at Cell Line Genetics® (WI), expression of Oct-3/4, SSEA-4, and SSEA-1 was evaluated using flow cytometry (BD FACSCalibur™) and directed differentiation into three germ layers was performed using human pluripotent stem cell functional identification kit (R&D Systems); nuclei were counterstained with Hoechst 33342 (blue).

Results


Cell morphology: Using NutriStem, cells displayed typical hPSC morphology on both surfaces.

Undifferentiated marker expression: hiPSCs expressed undifferentiation markers (Oct-3/4 and SSEA-4); differentiation marker (SSEA-1) was not detected.


Maintenance of pluripotency: hiPSCs were able to differentiate into three germ layers.

Karyotype: hiPSCs retained a normal karyotype.



Clump Passaging on Corning Matrigel Matrix (10 passages)

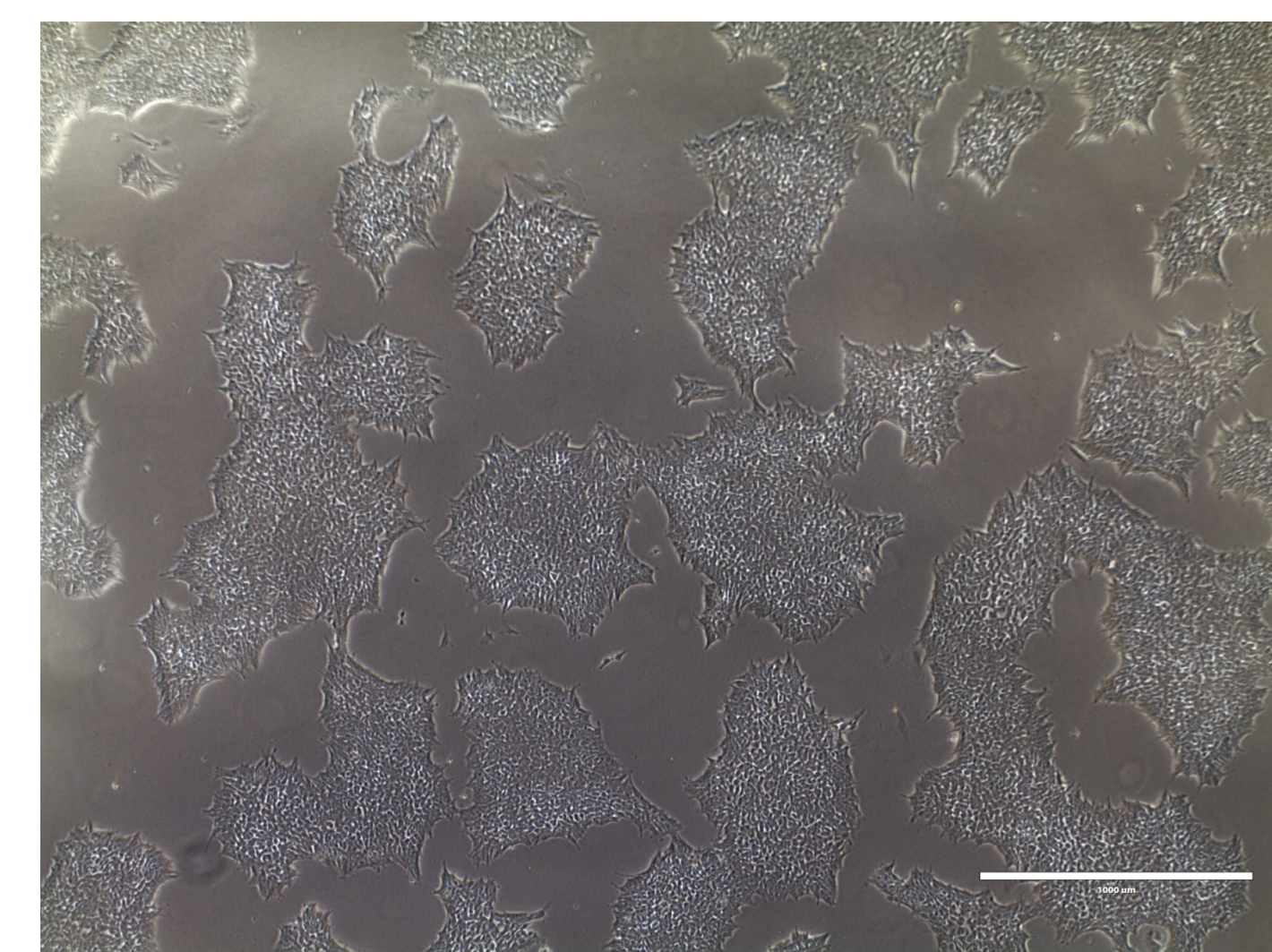


Morphology

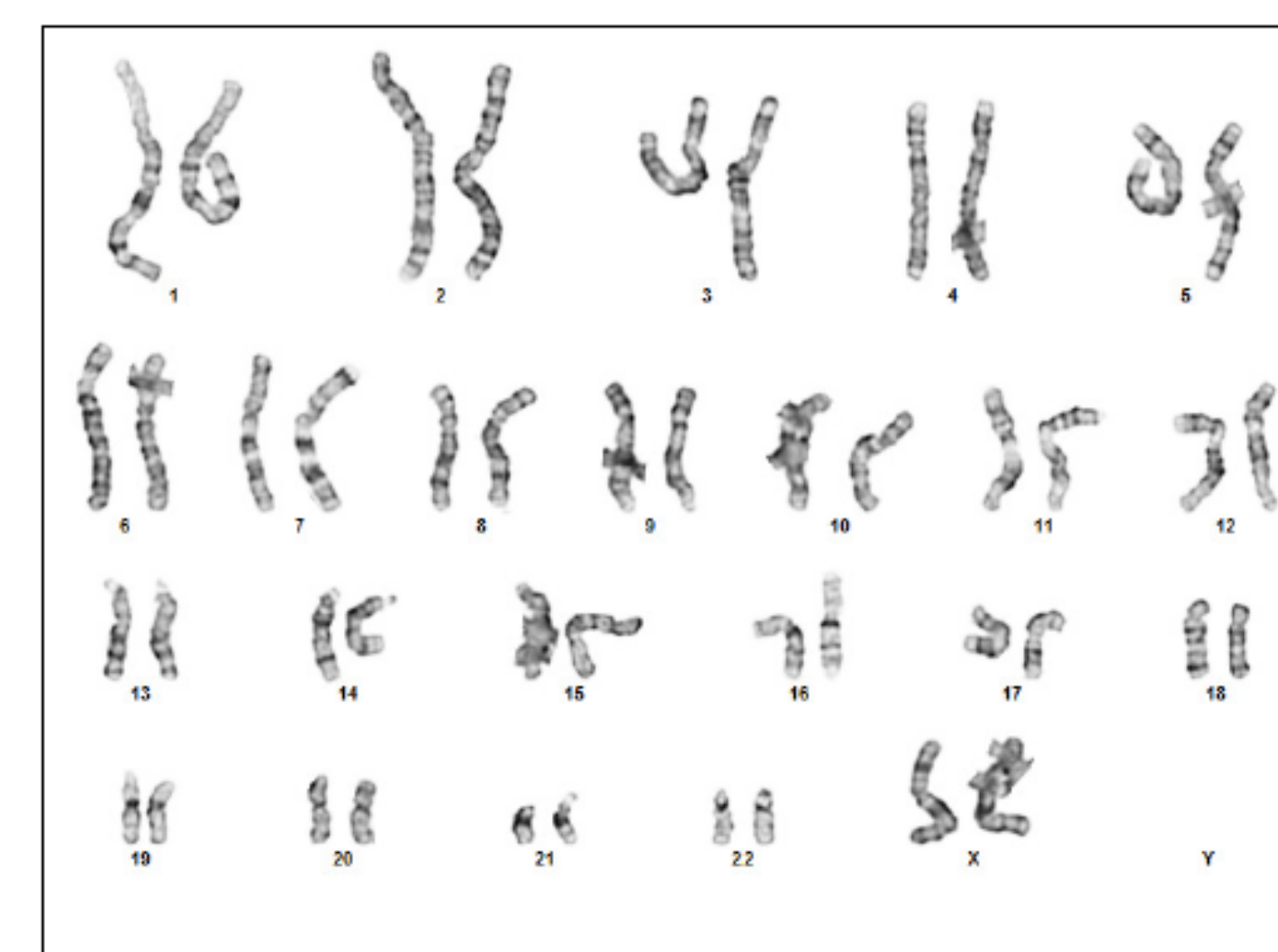


Karyotype

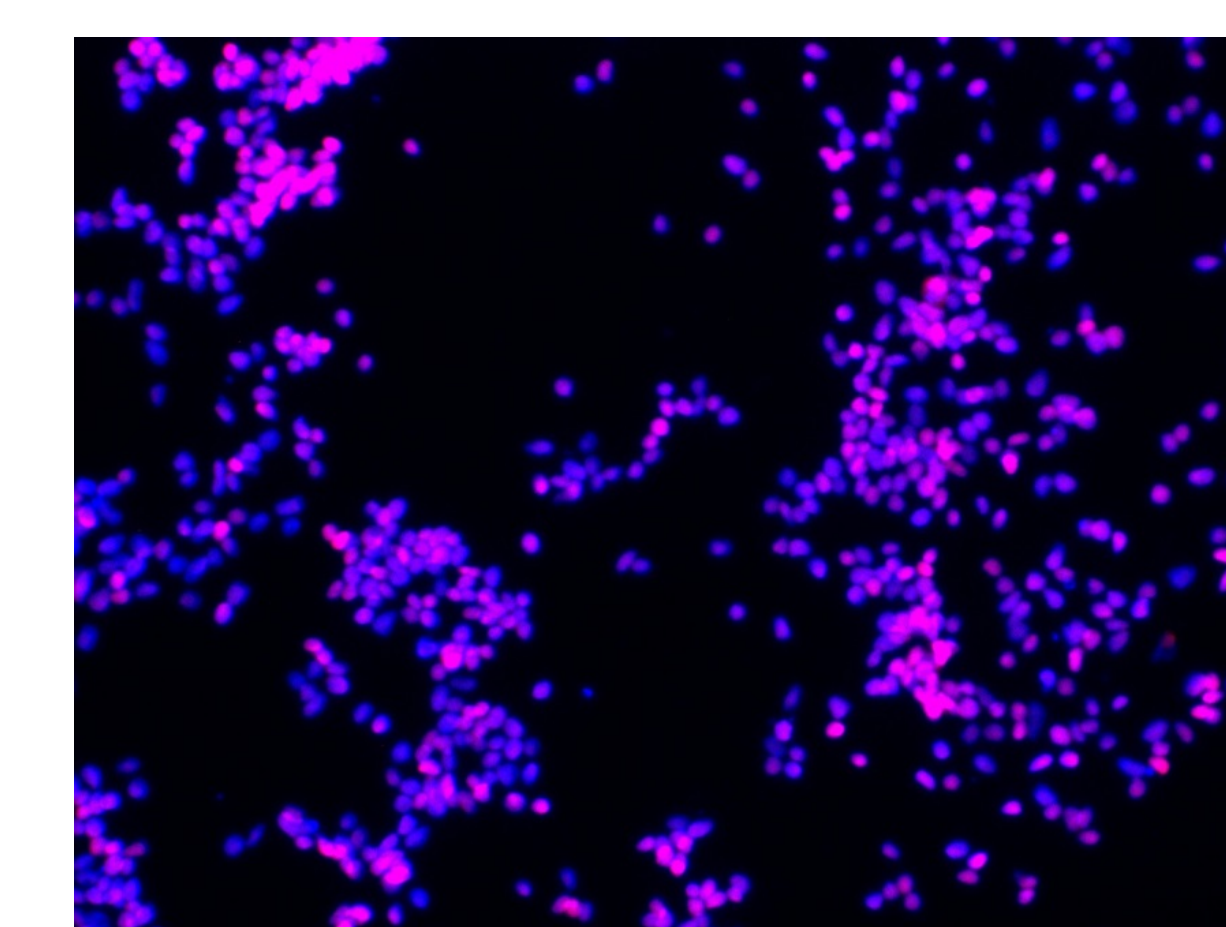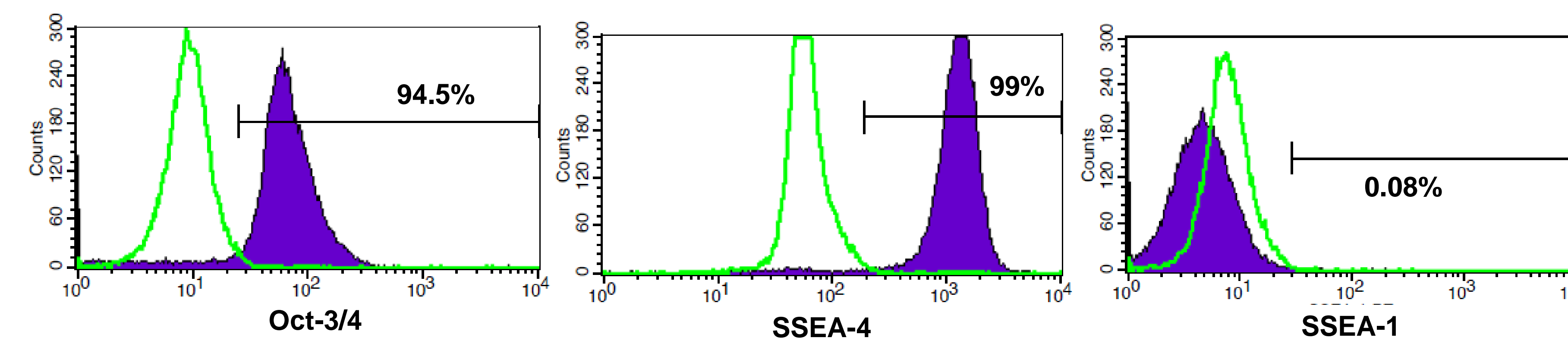

Endoderm (SOX17)

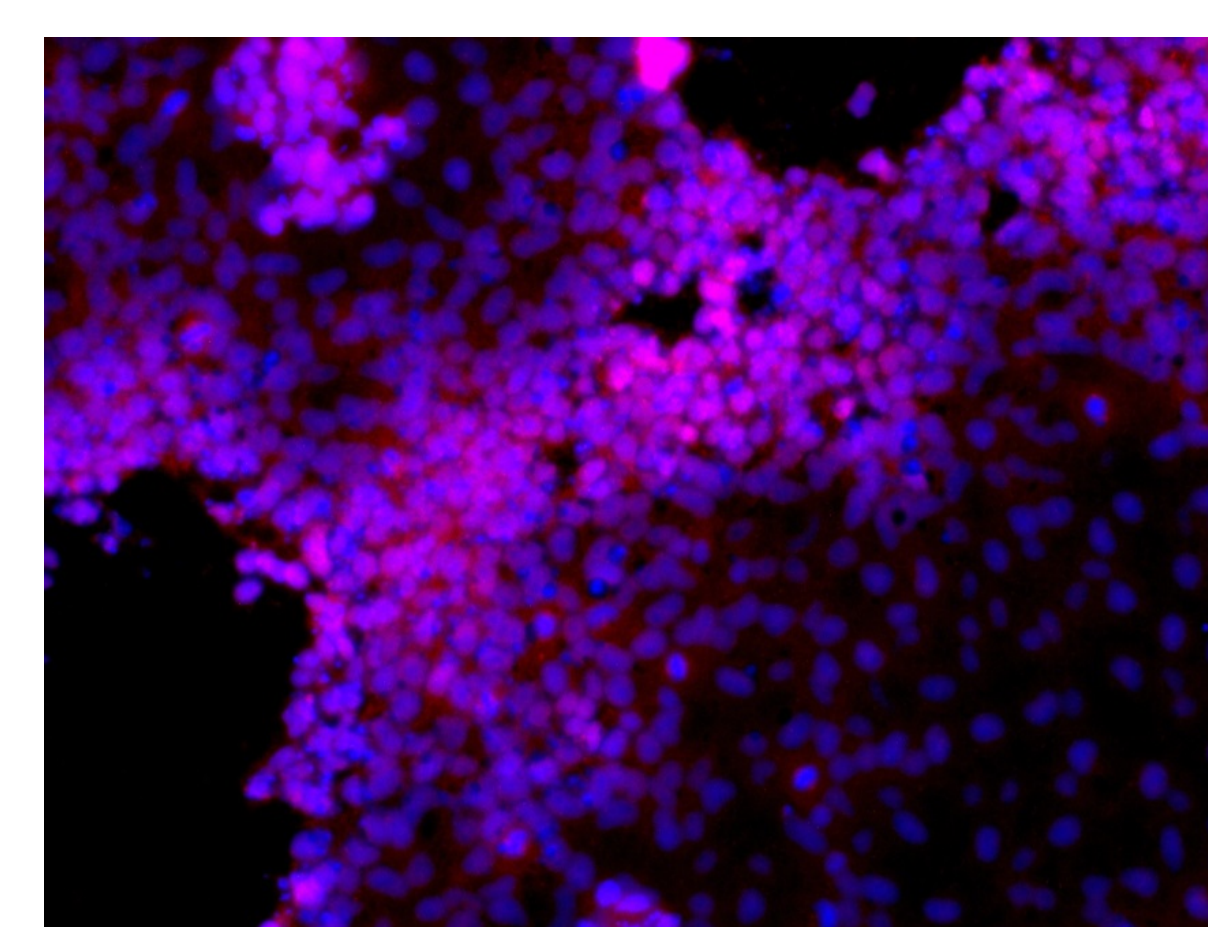


Ectoderm (Otx2)

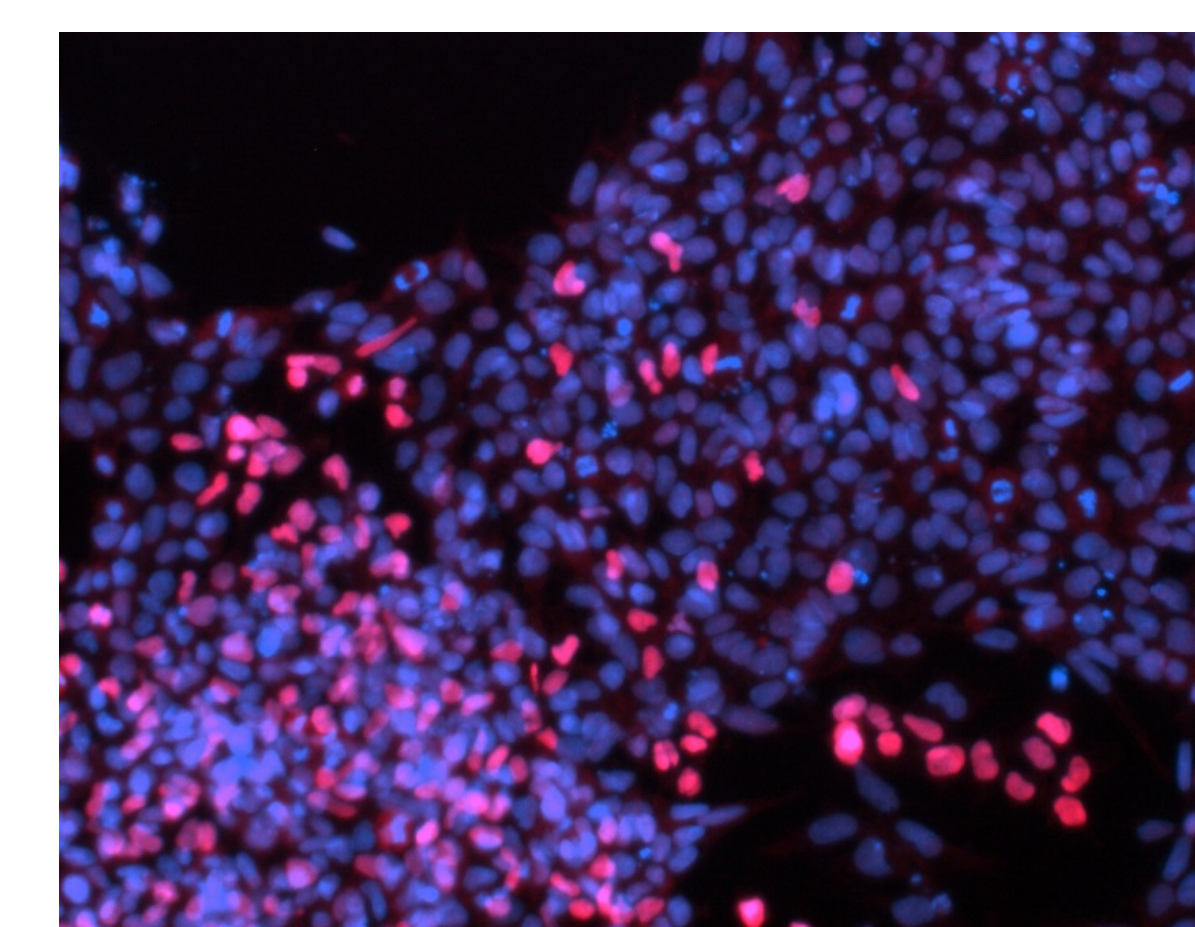


Mesoderm (Brachyury)



Single Cell Passaging on Corning PureCoat rLaminin-521 (10 passages)

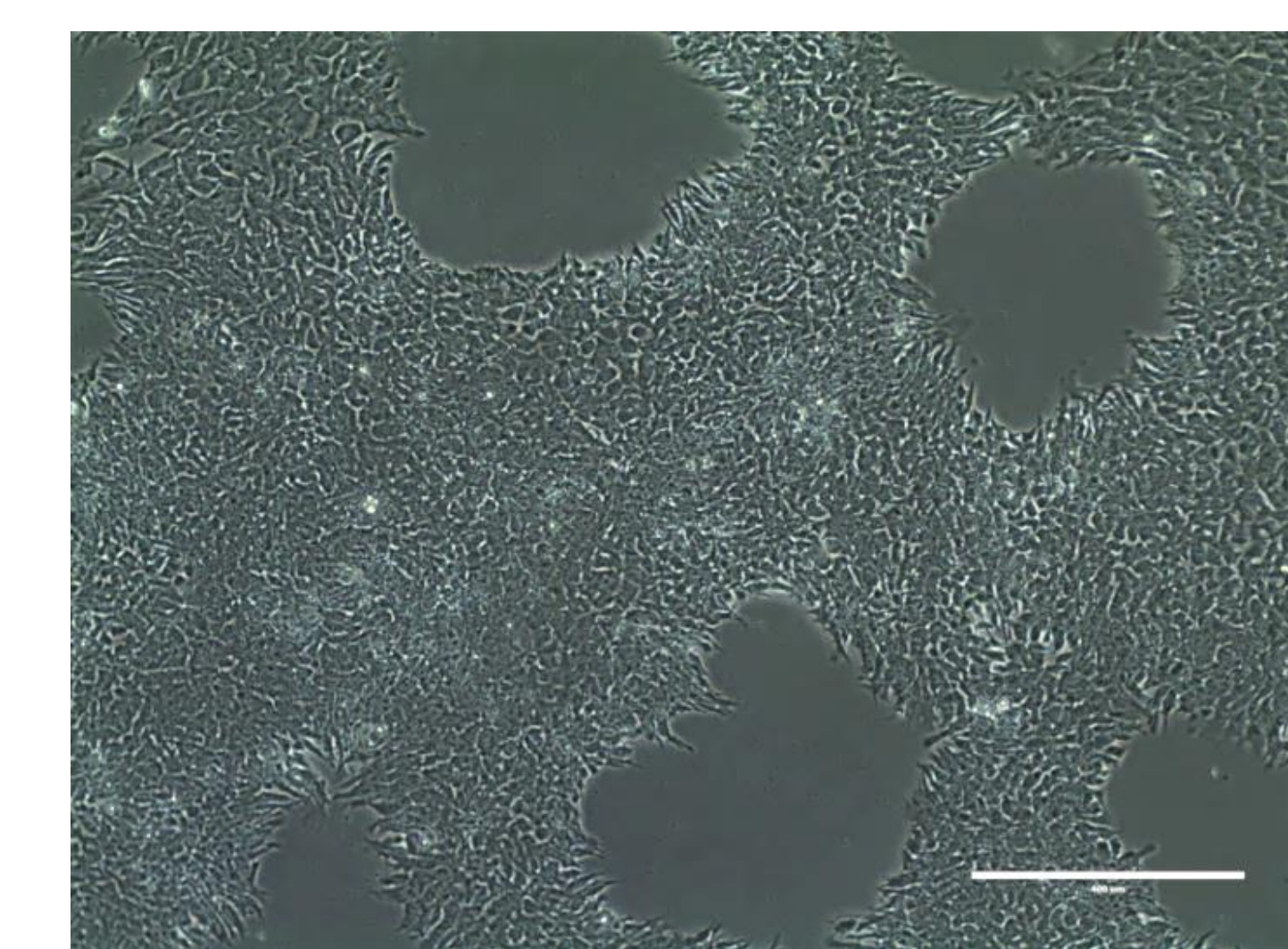


Morphology

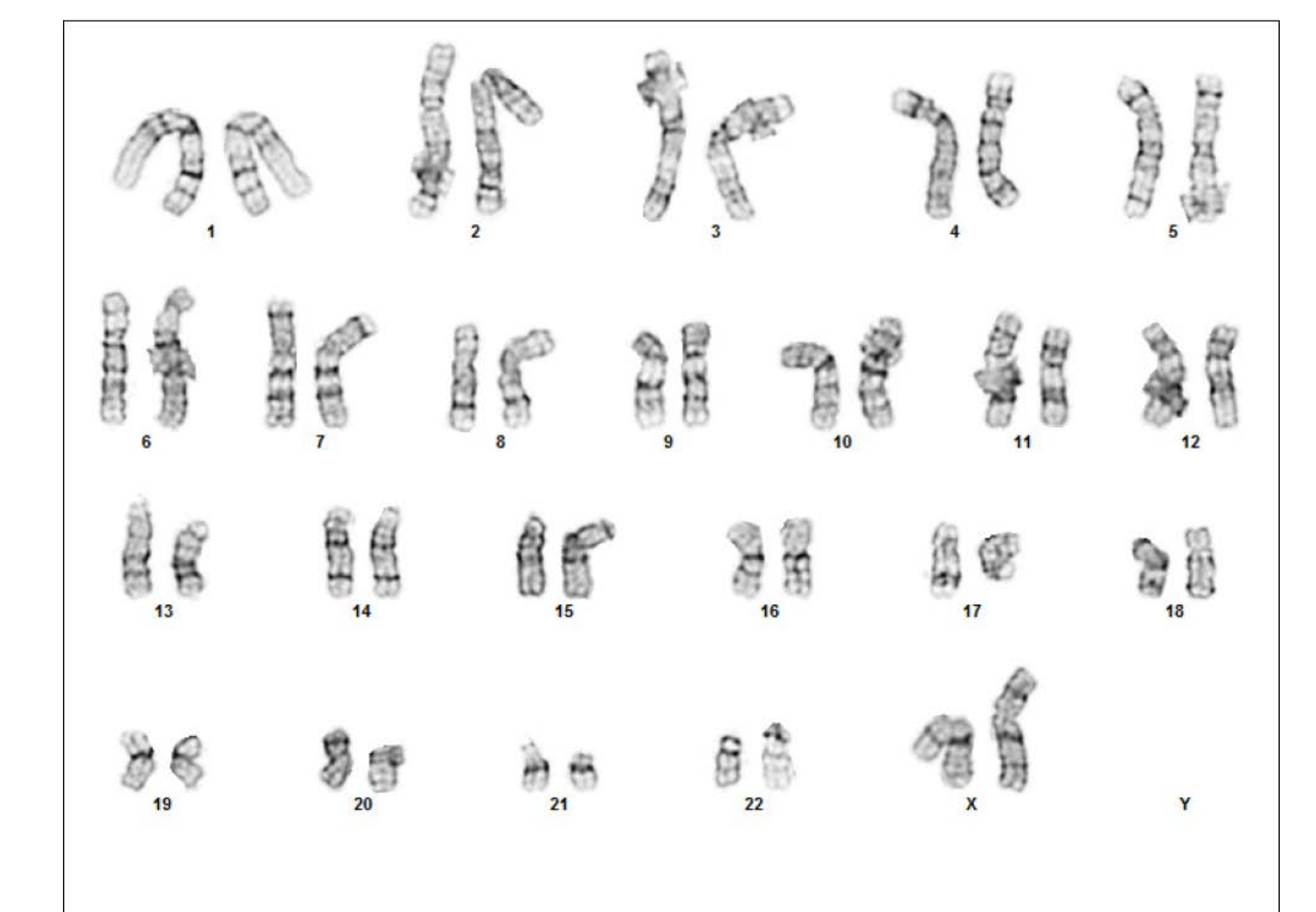


Karyotype

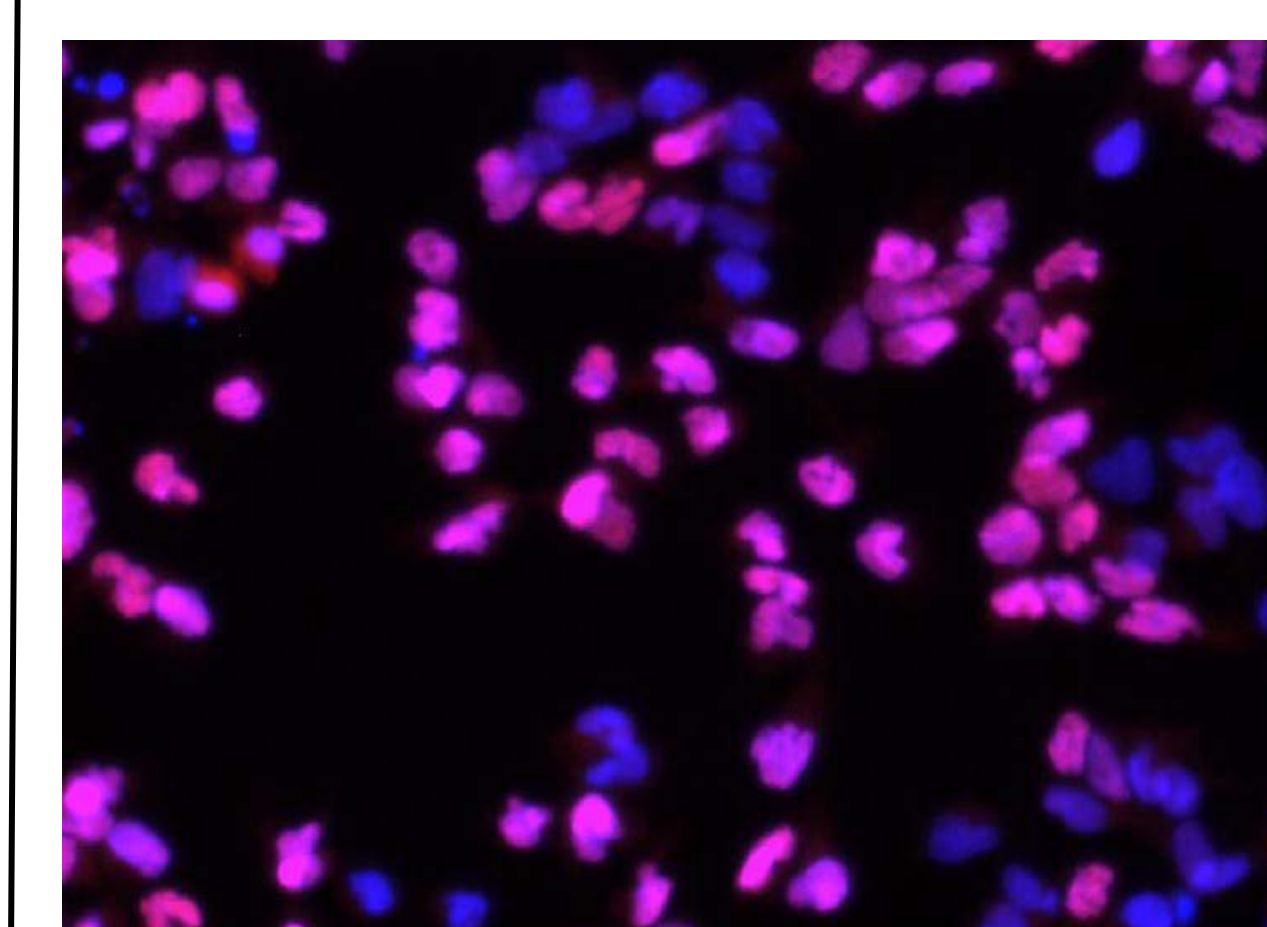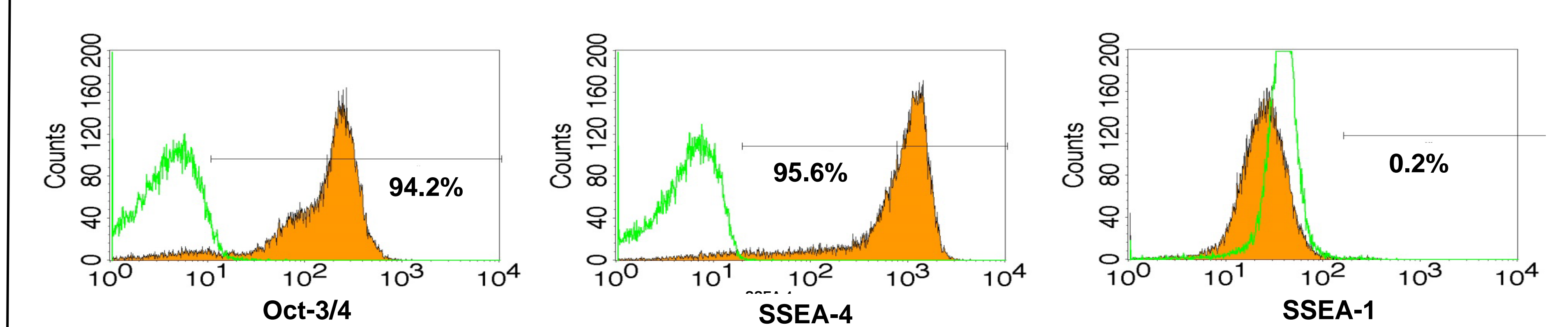

Endoderm (SOX17)

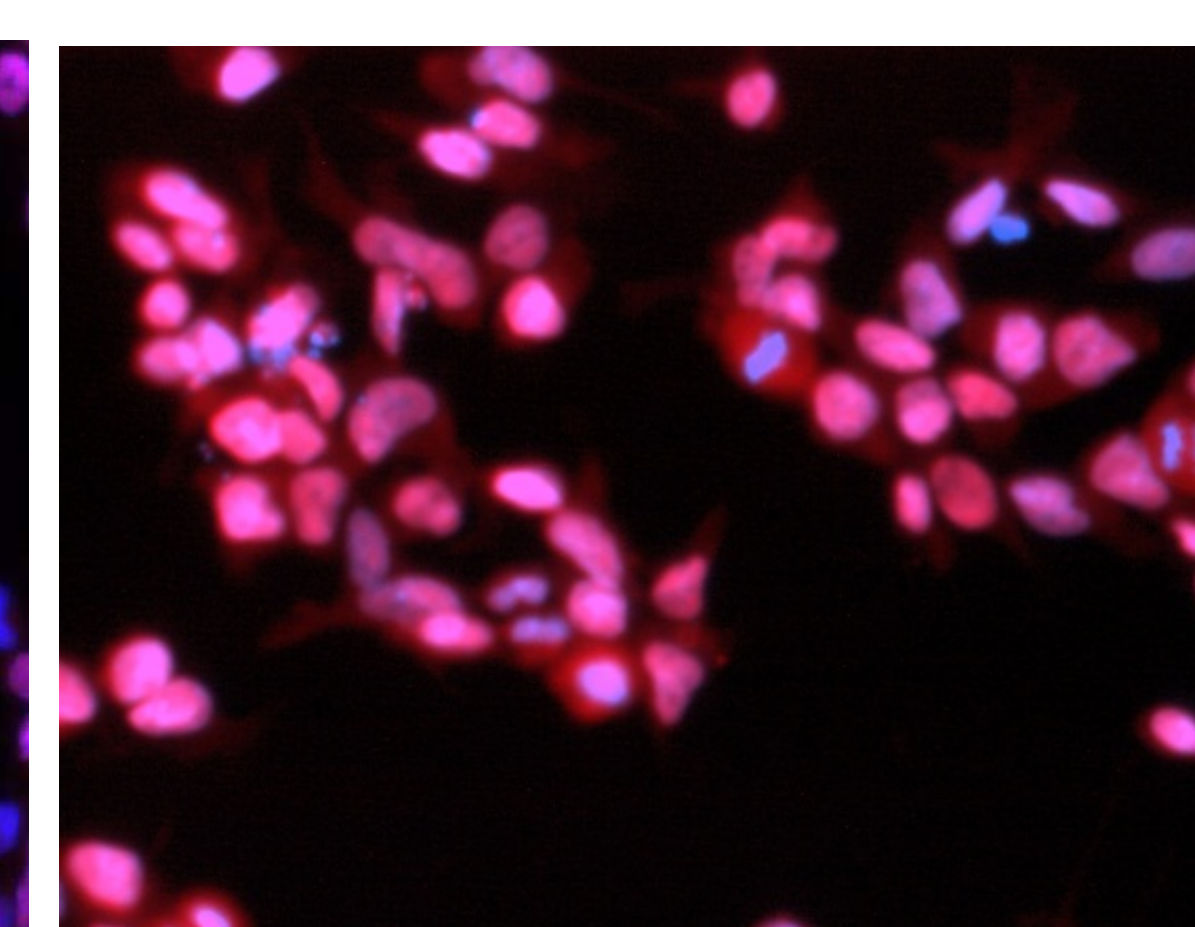

Ectoderm (Otx2)

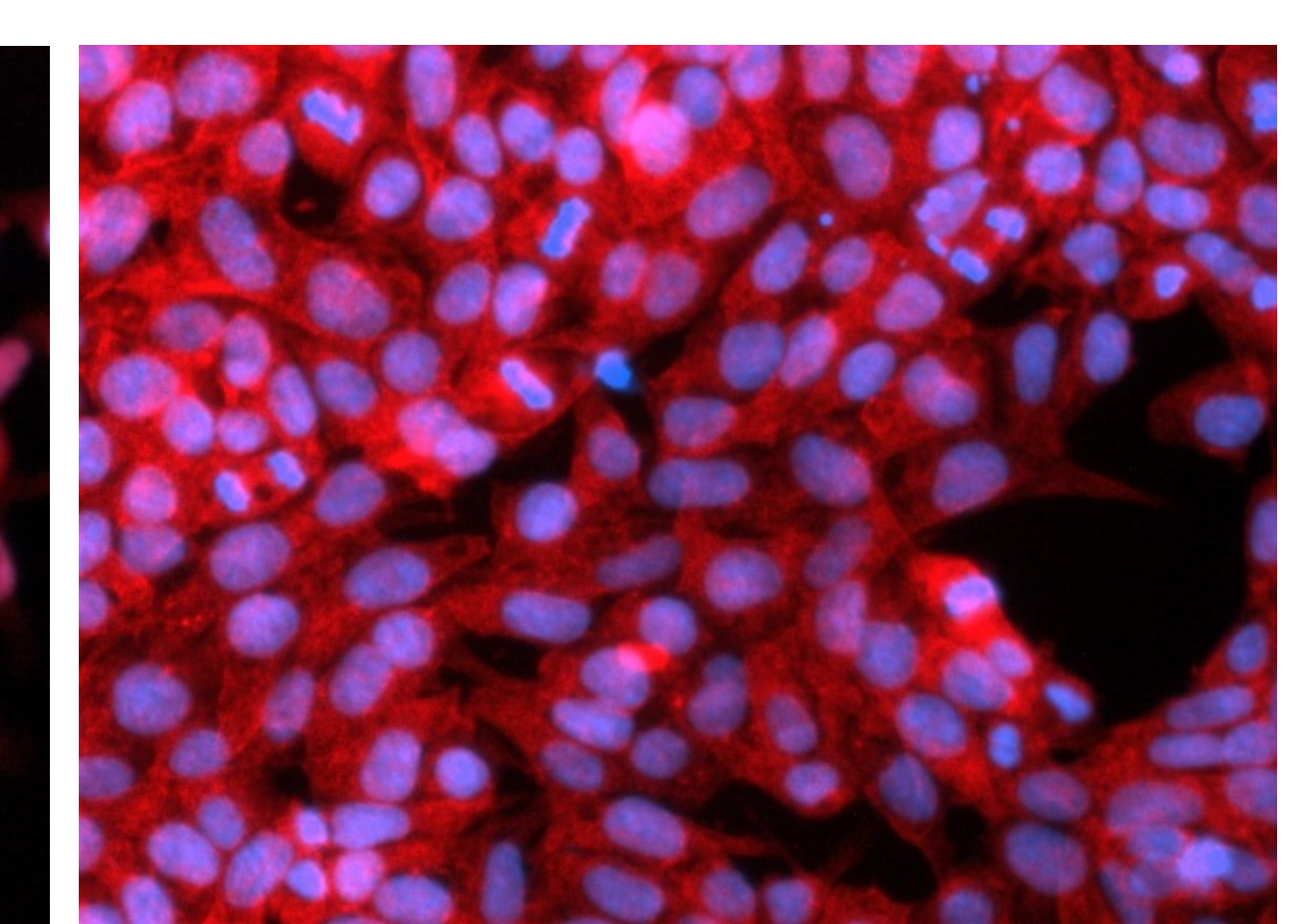

Mesoderm (Brachyury)

Conclusions

The xeno-free medium NutriStem allows usage of multiple surfaces and passaging methodologies for long-term culture and expansion of hPSCs.

- Clump passaging on Matrigel matrix
- Single cell passaging on PureCoat rLaminin-521 cultureware